

Technical Solutions

Optimisation of Spectrograph Dispersion Coefficients

Applicable for:

Andor Solis Software (version 4.24 onwards)
Shamrock Spectrograph Range


Andor Solis 2.24 introduces a new feature allowing the user to recalibrate the dispersion parameters following a change in the system (eg, changing camera or turret). The Dispersion Optimiser has been designed to fine-tune the dispersion coefficients ensuring more accurately labelled spectra on either side of the selected centre wavelength.

Using at least 2 known spectral peaks of a calibrated light source on each grating will ensure that the new values calculated are optimised for the useful wavelength ranges on all installed gratings. A good calibration will be obtained if 2 wavelengths on each grating are tracked, at a total of 6 if there are 3 gratings installed.


The wavelengths selected should be from either side of the wavelength region of interest, e.g. if a grating is being used to obtain spectra in the region of 400nm to 900nm then, using a Hg light source for example, the 912nm line should be used to ensure accuracy at the high end of the range and then either the 365nm line or the 435nm line in addition to cover the low end.

Note: This change is not reversible and therefore Andor recommend saving a copy of the original EEPROM before starting this process

The Dispersion Optimiser is found under the *Calibrate* menu in Solis.

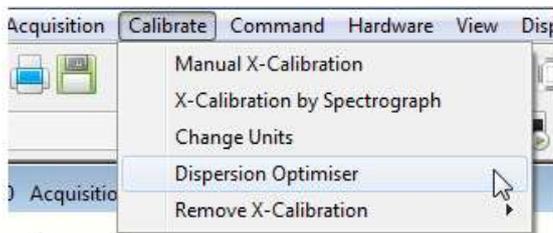


The Dispersion Optimiser menu is as follows:

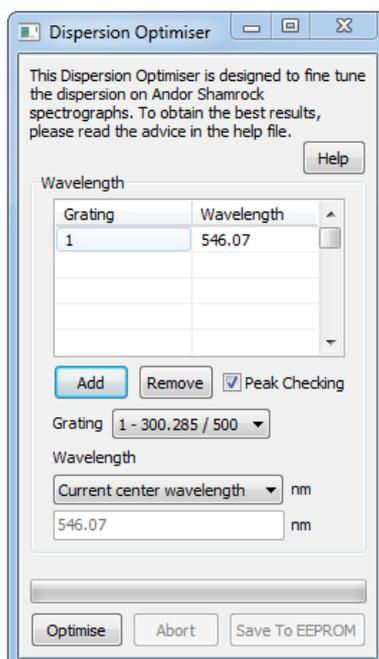


Steps to use the Dispersion Optimiser:

1) Select a Grating and Wavelength and add them to the list


- Select 1 grating, or the All Gratings Option:

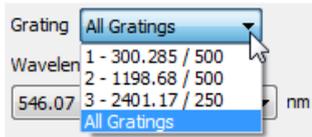

- Select a suitable wavelength:

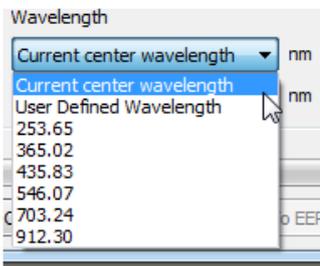


- Add the selected gratings and wavelengths to the list

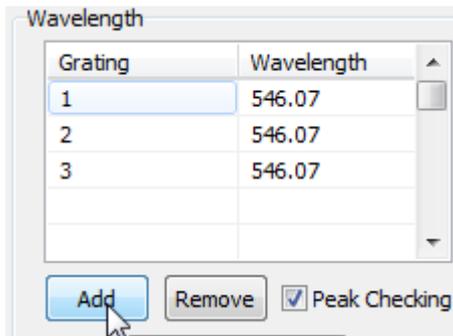


- Repeat for all gratings and wavelengths required

2) Click the *Optimise* button

- If *Peak Checking* is enabled, the following dialog will pop up at each of your chosen wavelength points to confirm the correct peak is being displayed

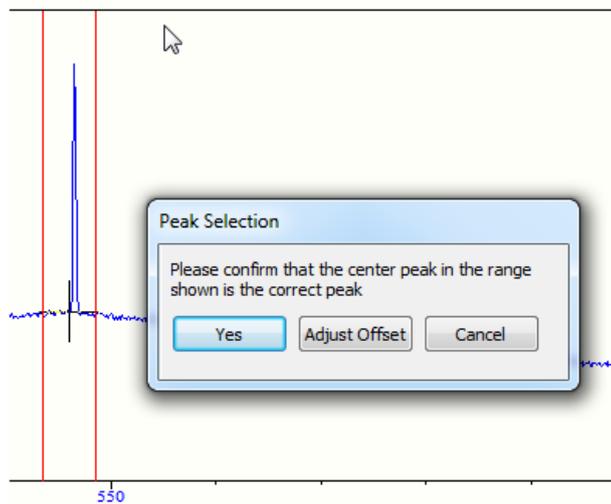


- Two red lines will be drawn around the centre range being used - the user can confirm that the peak they are expecting is within this range
- If the user selects 'Yes' the optimisation will continue
- If not in range, select *Adjust Offset*

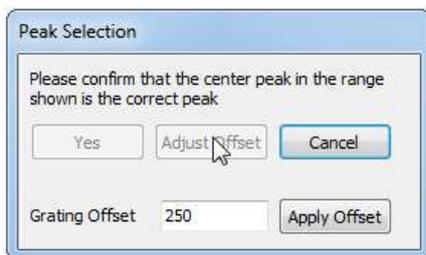

- By adjusting the grating offset, move the peak and until it is in range
- Selecting *Cancel* at any time will abort the current optimisation procedure

3) If satisfied with the new values, select *Save to EEPROM*

- Once the operation is complete the following display will be added to the dialogue

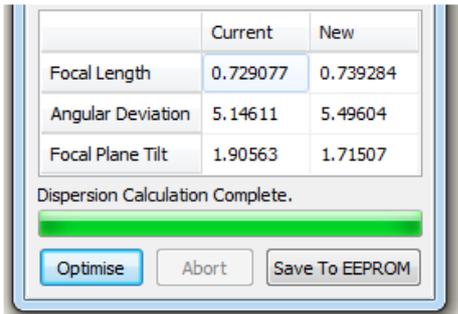


	Current	New
Focal Length	0.729077	0.739284
Angular Deviation	5.14611	5.49604
Focal Plane Tilt	1.90563	1.71507

Dispersion Calculation Complete.

Optimise Abort Save To EEPROM

- The 'Current' column shows the Optical Parameters currently in use, and the 'New' column shows the recommended values calculated by the Optimiser
- Select *Save to EEPROM* to update the new values to the EEPROM of the shamrock
- Once saving is complete the table of values will update to reflect the change in the current values

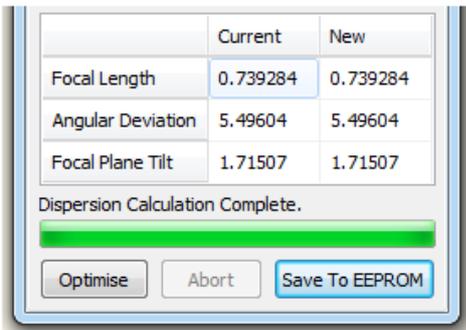

	Current	New
Focal Length	0.739284	0.739284
Angular Deviation	5.49604	5.49604
Focal Plane Tilt	1.71507	1.71507

Dispersion Calculation Complete.

Optimise Abort Save To EEPROM

For further assistance, please contact your local Andor Support Representative.