
iKon-M 934 Series

1 Megapixel, -100°C, 5 MHz Imaging CCD





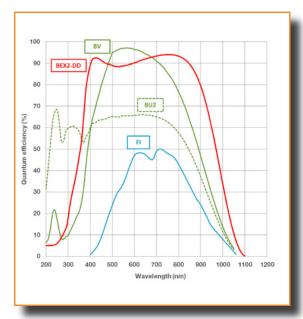

Features and Benefits

- Extended range Dual-AR option
 Superior UV-NIR broadband back-illuminated QE
- TE cooling to -100°C Critical for elimination of dark current required for long exposure experiments such as in vivo luminescence
- Fringe Suppression Technology™ NIR etaloning greatly reduced (BEX2-DD sensor types)
- Up to 5 MHz pixel readout ⁻⁸ High frame rates achievable
- Ultra low noise readout
 Intelligent low-noise electronics offer the most
 'silent' system noise performance available
- UltraVac^{™·1} Critical for sustained vacuum integrity and to maintain unequalled cooling and QE performance, year after year
- 13 x 13 μm pixel size Optimal balance of dynamic range and resolution
- Integrated shutter
 C-mount shutter as standard. Closed during
 readout to avoid vertical smear
- Cropped Sensor Mode
 Specialised acquisition mode for continuous
 imaging with fastest possible temporal resolution
- Enhanced Baseline Clamp
 Quantitative accuracy of dynamic measurements
- USB 2.0 connection
 Simple Plug & Play connection
- Integrated in EPICS
 Platform is fully integrated into the EPICS
 control software

Industry-Leading Ultra-Sensitive Imaging Technology

Andor's iKon-M 934 series cameras are designed to offer the ultimate in high-sensitivity, low noise performance, ideal for demanding imaging applications. These high resolution 1024 x 1024 CCD cameras boast up to 95% QE_{max} , high dynamic range, 13 µm pixels and exceptionally low readout noise. The iKon-M benefits from negligible dark current with industry-leading thermoelectric cooling down to -100°C. High sensitivity and the exceptionally low dark current makes iKon-M a perfect solution for the most demanding long exposure luminescence experiments such as plant imaging studies.

The 'Deep Depletion' sensor option offers ultimate performance for NIR applications, the new Dual AR coating extends the QE performance significantly across the UV/visible region of the spectrum for the broadest possible spectral coverage from one sensor. Fringe Suppression Technology™ radically minimizes etaloning effects in the NIR. Rapid vertical shifts combined with fast kinetics acquisition mode, comprehensive trigger modes and custom coated wedge window options, render the deep Depletion models ideal for NIR optimized Bose Einstein Condensation applications.

Specifications Summary

Active pixels	1024 x 1024
Sensor size	13.3 x 13.3 mm
Pixel size (W x H)	13 μm x 13 μm
Active area pixel well depth	100,000 e ⁻ (130,000 e ⁻ for BEX2-DD models)
Pixel readout rates (MHz)	5, 3, 1, 0.05
Read noise	2.9 e ⁻
Maximum cooling	-100°C
Frame rate	4.4 fps (full frame)

.



System Specifications²

Model number	DU934P	DU934P (Deep Depletion)
Sensor options	BU2: Back Illuminated CCD, UV-Enhanced, 250 nm optimized BV: Back Illuminated CCD, Vis optimized FI: Front Illuminated CCD	BEX2-DD : Back Illuminated CCD,Deep Depletion with fringe suppression, extended range dual AR coating
Active pixels	1024 :	x 1024
Pixel size	13 x 1	13 µm
Image area	13.3 x 13.3 mm w	ith 100% fill factor
Minimum temperatures ⁻³ Air cooled Coolant recirculator Coolant chiller, coolant @ 10°C, 0.75 l/min	-80°C -95°C -100°C	
Digitization	16-bit	
Blemish specifications	Grade 1 sensor from supplier. Camera blemishes as defined by Andor Grade A: www.andor.com/learning-academy/ccd-blemishes-and-non-uniformities-black-pixels-and-hot-pixels-on-a-ccd-sensor	
System window type	BU2 sensor: UV-grade fused silica, 'Broadband VUV-NIR', unwedged BV sensor: UV-grade fused silica, 'Broadband VIS-NIR', wedged BEX2-DD, FI sensors: UV-grade fused silica, 'Broadband VUV-NIR', wedged (other options available e.g. Bose-Einstein 780nm for Rb BEC experiments)	
Interface	USB 2.0	
Lens mount	C-mount	

Advanced Performance Specifications²

Dark current, e /pixel/sec *4 @ -80°C @ -100°C	0.00030 0.00012	0.01700 0.00047	
Pixel readout rates	5, 3, 1, 0.05 MHz		
Active area pixel well depth	100,000 e ⁻	130,000 e ⁻	
Read noise (e [.]) ⁻⁵ 0.05 MHz 1 MHz 3 MHz 5 MHz ⁻⁸	2.9 6.6 11.6 18.0	3.3 6.2 9.2 13.6	
Linearity *6	Better than 99%		
Vertical clock speed	11.3 to 67.3 μ s (software selectable)	4.25 to 64.25 μs (software selectable)	

Frame Rates⁷

		50 kHz		
Binning	Full Frame	512 x 512	256 x 256	128 x 128
1 x 1	0.04	0.1	0.2	0.4
2 x 2	0.2	0.2	0.4	0.8
4 x 4	0.6	0.6	0.9	1.5
8 x 8	2.0	1.2	1.7	2.8
16 x 16	5.1	2.3	3.2	5

		3 MHz		
Binning	Full Frame	512 x 512	256 x 256	128 x 128
1 x 1	2.6	5.2	10.0	18.6
2 x 2	6.3	10.9	19.3	33.2
4 x 4	13.4	20.9	34.2	53.5
8 x 8	25.1	36.3	54.4	76.5
16 x 16	41.5	56.1	76.5	97.1




5 MHz Visualization mode ⁻⁸				
Binning	Full Frame	512 x 512	256 x 256	128 x 128
1 x 1	4.4	8.5	16.0	28.8
2 x 2	8.4	15.9	28.5	47.6
4 x 4	15.6	28.1	47.0	70.8
8 x 8	27.3	45.8	69.4	93.5
16 x 16	43.6	66.9	91.2	111.5


1 Megapixel, -100°C, 5 MHz Imaging CCD



Quantum Efficiency Curves *

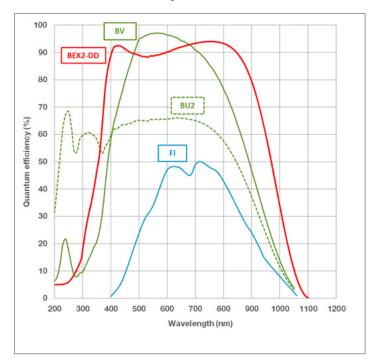

Application Image

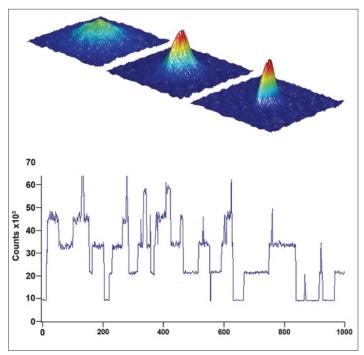


The top image shows surface plots of the transition from a thermal gas (left) to a Bose-Einstein condensate. The sharp, bimodal peak in the right figure is a signature of BEC. The graph below it shows fluorescence from a few-atom MOT against time, showing the discrete steps characteristic of single atoms entering and leaving the trap. Courtesy of Prof. Michael Chapman, Georgia Institute of Technology, USA.

Bioluminescence

High sensitivity and the exceptionally low dark current makes iKon-M a perfect solution for the most demanding long exposure luminescence experiments such as plant imaging studies. Read our full case study here.

Image Right: Example of bioluminescence timelapse of Arabidopsis to investigate the circadian clock. Courtesy of Professor Liu (Liu Lab, Shanghai Institutes for Biological Sciences, CAS, Shanghai)

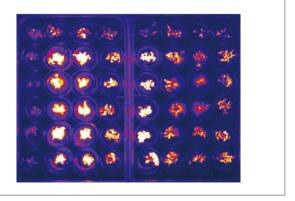

Have you found what you are looking for?

Need a larger sensor? The iKon-L 936 houses a 4 megapixel, back-illuminated CCD sensor (27.6 mm x 27.6 mm), cooling to -100°C, low noise performance and up to 5 MHz readout.

Need the ultimate in sensitivity? The iXon back-illuminated EMCCD series offers > 90% QE and single photon sensitivity, combined with fast frame rate performance.

Need faster frame rates? The Neo and Zyla sCMOS deliver up to 100 frames/sec (full frame).

Need smaller pixels? Check out the the Neo and Zyla sCMOS.

Need a customised version? Please contact us to discuss our Customer Special Request options.

Check out Andor's New Sona sCMOS. back-illuminated sCMOS delivers ultra-sensitivity, high speed, high-resolution, large field of view & high dynamic range!

3



Creating The Optimum Product for You





Step 1.	Choose the sensor type option	
	Description	Code
3 3	Back Illuminated CCD, Vis optimized	BV
	Front Illuminated CCD	FI
Sensor	Back Illuminated CCD, AR coated for optimized performance in the 250 nm region	BU2
Туре	Back Illuminated CCD, Deep Depletion with fringe suppression, extended range dual AR coating	BEX2-DD

Step 2. Select an alternative camera window (optional)

The standard window has been selected to satisfy most applications. However, other options are available. The alternative camera window code must be specified at time of ordering.

To view and select other window options please refer to the 'Camera Windows Supplementary Specification Sheet' which gives the transmission characteristics, product codes and procedure for entering the order. Further detailed information on the windows can be found in the Technical note - 'Camera Windows: Optimizing for Different Spectral Regions'.

Step 3. Select the required accessories and adapters Description Order Code XW-RECR Re-circulator for enhanced cooling performance (supplied with 2x2.5 m tubing as standard) Oasis 160 Ultra compact chiller unit (tubing to be ordered separately) ACC-XW-CHIL-160 ACC-6MM-TUBING-2X2.5/ 6 mm tubing options for ACC-XW-CHIL-160 (2x2.5 m or 2x5m lengths) ACC-6MM-TUBING-2X5M C-mount to Canon FD-mount adapter OA-CCFM OA-CNAF C-mount to Nikon F-mount adapter Accessories & Adapters C-mount to Olympus OM adapter OA-COFM Auto extension tubes (set of 3) for C-mount OA-ECMT Auto extension tubes (set of 3) for Nikon F **OA-ENAF** ACC-USBX-EU USB Extender: Icron USB 2.0 Ranger (supports up to 100 m) - EU/UK/US ACC-USBX-UK ACC-USBX-US 30 m Ethernet cable (for use with the above ACC-USBX-** USB extenders) ACC-ELC-13295

Step 4. Select the required software

display and export.

The iKon-M requires at least one of the following software options:

Software

Camera

Window

Andor SDK A software development kit that allows you to control the Andor range of cameras from your own application. Available as 32/64-bit libraries for Windows (8, 8.1 and 10), compatible with C/C++, C#, Delphi, VB.NET, LabVIEW and Matlab. Linux SDK compatible with C/C++.


Third party software compatibility Drivers are available so that the iKon-M range can be operated through a large variety of third party imaging packages. See Andor web site for detail: https://andor.oxinst.com/learning/view/article/third-partyimaging-software-support

Solis for Imaging A 32-bit and fully 64-bit enabled application for Windows (8, 8.1 and 10) Linux and Labview, offering rich functionality for data acquisition and processing. AndorBasic provides macro language control of data acquisition, processing,



Product Drawings

Dimensions in mm [inches]

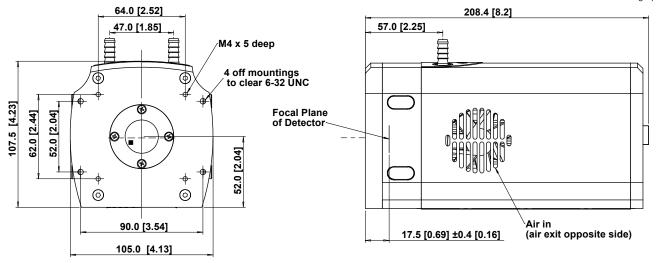

■ = position of pixel 1,1 Weight: 2.2 kg [4 lb 13 oz]

Connecting to the iKon-M

Camera Control Connector type: USB 2.0 with optional lockable connector

TTL / Logic

Connector type: SMB, provided with SMB - BNC cable Fire (Output), External Trigger (Input), Shutter (Output)


I²C connector

Compatible with Fischer SC102A054-130 Shutter (TTL), I²C Clock, I²C Data, +5 Vdc, Ground


Minimum cable clearance required at rear of camera 90 mm

Typical Applications

- Astronomy
- Bioluminescence/Chemiluminescence
- **Biochip Reading**
- Bose-Einstein Condensation (BEC)
- Photovoltaic Inspection (see the iKon-M PV Inspector Specification Sheet)
- Semiconductor analysis

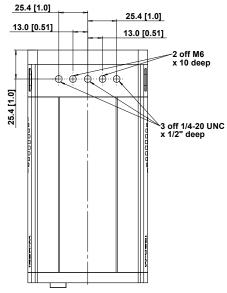

Mounting hole locations

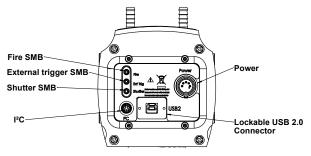

Rear connector panel

Third

iKon-M 934 Series

1 Megapixel, -100°C, 5 MHz Imaging <u>CCD</u>





Order Today

Need more information? At Andor we are committed to finding the correct solution for you. With a dedicated team of technical advisors, we are able to offer you one-to-one guidance and technical support on all Andor products. For a full listing of our regional sales offices, please see: www.andor.oxinst.com/contact

Our regional headquarters are:

Europe Belfast, Northern Ireland Phone +44 (28) 9023 7126 Fax +44 (28) 9031 0792

North America Concord, MA, USA Phone +1 (860) 290 9211 Fax +1 (860) 290 9566 Japan Tokyo Phone +81 (3) 6732 8968 Fax +81 (3) 6732 8939

China Beijing Phone +86 (10) 5884 7900 Fax +86 (10) 5884 7901

Footnotes: Specifications are subject to change without notice

- Assembled in a state-of-the-art cleanroom facility, Andor's UltraVac[™] vacuum process combines a permanent hermetic vacuum seal (no o-rings), with a stringent protocol to minimize outgassing, including use of proprietary material.
- 2. Figures are typical unless otherwise stated.
- Specified minimum air cooled temperature assumes ambient temperature of 25°C. Specified minimum temperature with coolant assumes coolant temperature of 10°C.
- 4. The dark current measurement is averaged over the sensor area excluding any regions of blemishes.
- Readout noise is for the entire system. It is a combination of sensor readout noise and A/D noise. Measurement is for Single Pixel readout with the sensor at a temperature of -80°C and minimum exposure time under dark conditions.
- 6. Linearity is measured from a plot of counts vs exposure time under constant photon flux up to the saturation point of the system.
- The frame rates shown are for the BEX2-DD model, for a range of binning or array size combinations. All
 measurements are made with 4.25 µs vertical shift speed. It also assumes internal trigger mode of operation
 and minimum exposure time.
- 8. 5 MHz is for focusing/visualization mode only.
- 9. Quantum efficiency of the sensor at 25°C as supplied by the sensor manufacturer.



Minimum Computer Requirements:

Items shipped with your camera

1x 2 m BNC - SMB connection cable

1x Power supply with mains cable

1x CD containing Andor user guides

1x 3 m USB 2.0 cable Type A to Type B

1x Individual system performance booklet

- 3.0 GHz single core or 2.4 GHz multi core processor
 2 GB BAM
- 100 MB free hard disc to install software (at least 1GB recommended for data spooling)
- USB 2.0 High Speed Host Controller capable of sustained rate of 40MB/s
- Windows (8.1 and 10) or Linux

Operating & Storage Conditions

- Operating Temperature: 0°C to 30°C ambient
- Relative Humidity: < 70% (non-condensing)
- Storage Temperature: -25°C to 50°C

Power Requirements

- 100 240 VAC, 50 60 Hz
- Power consumption: 48W max.



Windows is a registered trademark of Microsoft Corporation. LabVIEW is a registered trademark of National Instruments. MATLAB is a registered trademark of The MathWorks Inc.