

# iXon Ultra


# The World's Highest Performance Back-illuminated EMCCDs

## **Key Specifications**

- ✓ Single photon sensitive
- ✓ Exceptional QE > 95%
- ✓ 13 or 16 µm pixel size
- ✓ Active pixels: 1024 x 1024 or 512 x 512
- ✓ TE cooling down to -95 or -100°C
- ✓ 26 or 56 fps full frame
- ✓ SRRF-Stream<sup>+</sup> real time super-resolution

## **Key Applications**

- Quantum imaging
- ✓ Cold atom & ion research
- ✓ Fast astronomy
- ✓ Tomography
- √ Fast spectroscopy
- ✓ Single molecule detection
- ✓ Super-resolution

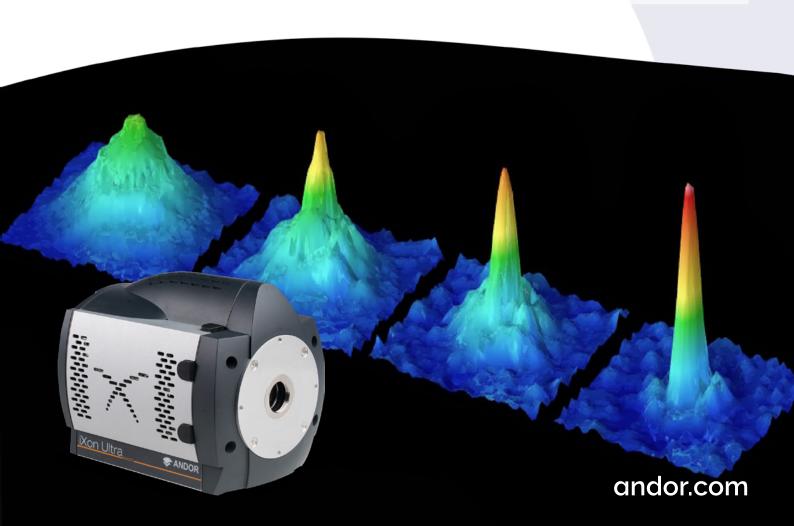

## iXon Ultra

## iXon Ultra 888: Field of View & Sensitivity... Now 3x Faster!

The highly innovative **iXon Ultra 888** megapixel, back-illuminated EMCCD camera offers single photon sensitivity across a large field of view, at 26 fps. Building on a rich history of first to market innovation, the 'supercharged' iXon Ultra 888, represents a massive performance boost for the largest available EMCCD sensor, as well as the first USB 3.0 enabled EMCCD camera.

The iXon Ultra 888 has been fundamentally re-engineered to facilitate a 3x acceleration of the pixel readout speed to an unprecedented 30 MHz, whilst maintaining quantitative stability, propelling the full frame performance to video rate. Furthermore, Andor's unique 'Crop Mode' can be employed to further boost frame rates from a user defined sub-region, for example pushing a 512 x 512 sub-array to 93 fps and a 128 x 128 area to 697 fps.

With a 1024 x 1024 sensor format and 13  $\mu$ m pixel size, the resolving power, field of view and unparalleled speed of the iXon Ultra 888 render it the most attractive and versatile EMCCD option for demanding applications such as single molecule detection, super-resolution microscopy, live cell imaging and high time resolution astronomy.

The iXon Ultra is designed to be the most flexible yet easy to use EMCCD on the market, optimizable for a wide variety of application requirements in a single click via the OptAcquire™ feature. Furthermore, signal can be quantitatively calibrated in units of electrons or photons, either in real time or post-processing. Patented, pioneering technology offers automated recalibration of EM gain, alongside anti-ageing protection.


Crucially, the iXon brand carries an outstanding reputation within the industry for quality and reliability, brandishing an unparalleled track record of minimal field failures.



Image comparison of a fluorescently labelled fixed BPAE cell, recorded with a widefield fluorescence microscope and a SRRF-Stream enabled iXon 888 camera.

## **Key Specifications**

Active pixels (H x V)	1024 x 1024
Pixel size (W x H; µm )	13 x 13
Image area (mm)	13.3 x 13.3
Active Area Pixel Well Depth (e-)	80,000
Max Readout Rate (MHz)	30
Frame rates (fps)	26 (full frame) - 9690
Read noise (e-)	<1 with EM gain
QE Max	>95%

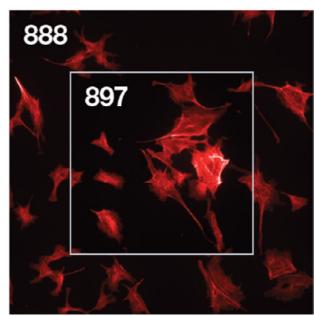

Field of View Comparison between iXon Ultra models. The 888 model has a  $\times 2.6$  greater sensitive area than the 897 model.

## iXon Ultra 897:

## The market leading backilluminated EMCCD, now accelerated to 56 fps.

The iXon Ultra 897 platform takes the popular back-illuminated 512 x 512 frame transfer sensor and overclocks readout to 17 MHz, pushing speed performance to an outstanding 56 fps (full frame), whilst maintaining single photon sensitivity and quantitative stability throughout. New Optically Centred Crop Mode unlocks unparalleled frame rate performance from centrally located ROIs, ideal for the particular speed and sensitivity requirements of super-resolution microscopy.

The iXon Ultra maintains all the advanced performance attributes that have defined the industry-leading iXon range, such as deep vacuum cooling to -100°C, extremely low spurious noise, and Andor's patented EM gain recalibration technology (EMCA™). Count Convert functionality allows real time data acquisition in units of electrons or incident photons and OptAcquire facilitates one-click optimization of this versatile camera to a variety of application conditions.


Additional features of the iXon Ultra include plug and play USB connectivity, a lower noise conventional CCD mode and an additional Camera Link output, offering the unique ability to directly access data for 'on the fly' processing, ideally suited to data intensive applications such as adaptive optics or superresolution microscopy.



## **Key Specifications**

Active pixels (H x V)	512 x 512
Pixel size (W x H; µm )	16 x 16
Image area (mm)	8.2 x 8.2
Active Area Pixel Well Depth (e-)	180,000
Max Readout Rate (MHz)	17
Frame rates (fps)	56 (full frame) - 11,074
Read noise (e-)	<1 with EM gain
QE Max	>95%

## Stability Plot

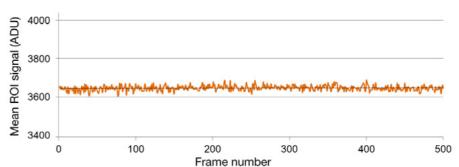

EM Gain stability in the iXon Ultra 897 @ 55 fps. 500 frame kinetic series; frame transfer (overlapped) acquisition; 17.8 ms exposure time; x300 EM gain.

---

-

## Features & Benefits

Single Photon Sensitive & > 95% QE	Optimal SNR in light starved applications such as single molecule detection and quantum physics.
NEW 'SRRF-Stream*' (optional)	Real time, cell super-resolution functionality. Living and fixed cells, works on most modern fluorescence microscopes. Super-resolution down to 50 nm.
Blue optimized model	Our new sensor option offers up to 20% higher sensitivity to photons across the blue region of the spectrum, ideal for quantum and ion imaging as well as blue-emitting BEC.
Overclocked readout speeds	Follow dynamic changing processes.
Crop Mode	Continuous imaging with fastest possible frame rate from centrally positioned ROIs. Highly enabling for live cell super-resolution and much more (e.g. 251 fps with 256 $\times$ 256 ROI).
TE cooling to -100°C	Elimination of dark current detection limit.
RealGain™	Absolute EMCCD gain selectable directly from a linear and quantitative scale.
Lower Noise CCD Mode	'2 in 1' flexibility. EMCCD for ultra-sensitivity at speed, conventional CCD for longer acquisitions.
Fringe Suppression (optional)	Reduced etaloning in NIR.
OptAcquire	Optimize the highly flexible iXon for different application requirements at the click of a button.
Count Convert	Quantitatively capture and view data in electrons or incident photons. Count Convert does this important conversion for you.
EMCAL™	Patented user-initiated self-recalibration of EM gain.
Qualified down to -20°C ambient temperature	Excellent for use at observatories.
Minimal Clock-Induced Charge	Confident discrimination of single photon events in Quantum Imaging.
UltraVac™	Critical for sustained vacuum integrity and to maintain unequalled cooling and QE performance, year after year. Seven year vacuum warranty.
Spurious Noise Filter	Intelligent algorithms to filter clock induced charge events from the background.
Direct Data Access	Camera Link output port to facilitate direct access to data for 'on the fly' processing.
Enhanced photon counting modes	Intuitive single photon counting modes ideal for Quantum Imaging. Real time or post-processing.
FPGA Timestamp	Hardware generated timestamp with 10 ns accuracy.
ASTRO.control compatibility (Ultra only)	Redlogix ASTRO.control is a dedicated platform for control of astronomical telescopes and instrumentation <a href="http://www.andor.com/astrocontrol.aspx">http://www.andor.com/astrocontrol.aspx</a>

# **Application Focus**

## The Physicist's Choice

The unique high-performance specifications of the optimized iXon range have been serving the physical scientist and astronomer in scenarios that demand more than simply an EM sensor in a camera. Andor have worked with numerous scientists to deliver solutions that work for their particular application requirements, such as specific coatings, coupling to fibre optic scintillators and 'minimized latency' direct data access solutions.

## Quantum Science

Quantum entanglement is now being studied with the aim of providing insight into the practical applications of this phenomenon. Quantum cryptography, communication and computing may soon rely on high-fidelity readouts of entangled photons. To detect these single particles researchers must rely on the most sensitive detection as is provided by iXon Ultra EMCCD cameras.

## Cold Atom & Ion Research

Our new blue-enhanced back-illuminated EMCCD sensor option is ideal for cold atom imaging experiments such as blue-emitting Bose-Einstein Condensation (BEC), ion traps and degenerate matter studies. Single photon sensitivity enables users to discern signals from single trapped ions and atoms.

## Fast Astronomy

Extremely fast and ultra-sensitive performance of binned sub-regions make the iXon Ultra an ideal Shack Hartmann AO detector. The large FOV of the Ultra 888 and fast frame rate are ideal for Lucky / Speckle Imaging. Qualified down to -20°C ambient, perfect for night observation.

## Tomography


Lens coupled or customer fibre coupled, the iXon Ultra provides the superlative ultrasensitive, large array solution for fast tomography.

## Fast Spectroscopy


The iXon Ultra has been a very valuable detector for fast spectroscopy applications, such for rapid spectroscopic mapping or fast monitoring of fast chemical reactions. When operated in full vertical binning the cameras can be pushed to thousands of spectra per second, accessing sub-ms dynamics. The iXon Ultra camera models are fully compatible with Andor's Kymera and Shamrock range of high-end spectrographs.

## iXon Ultra 888 - Capture everything

- Extreme Sensitivity capture, resolve and quantify extremely weak structures.
- ✓ **Capture More** the 13.3 x 13.3 mm sensor of the Ultra 888 is well suited to the optically useful FOV from a microscope.
- 3x Faster more and more studies of cell processes require greater temporal resolution. The Ultra 888 offers the fastest available EMCCD readout capability.
- Superb Image Quality megapixel resolution and uniform backgrounds for publication quality imaging.
- ✓ 2 in 1 the Ultra 888 can operate as an EMCCD or a low noise CCD, rendering it optimal for both fast and slow capture.

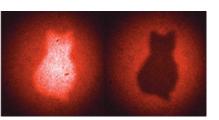


Quantum entanglement, see Quantum imaging finally saves Schrödinger's cat. Courtesy of Anton Zeilinger, Institute for Quantum Optics and Quantum Information, University of Vienna.

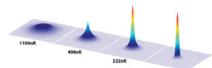

A phase transition and BEC formation in a gas of erbium atoms imaged at 401 nm matched by camera's QE at approx. 75%. Courtesy of K. Aikawa et al., University of Innsbruck, Austria.



The iXon Ultra 888 serves as the Focal Plane Imager of the SOFIA telescope. Courtesy of Pasquale Temi & E.E. Becklin, NASA.

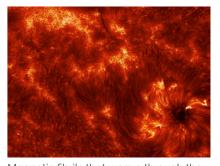

Magnetic fibrils that weave through the solar chromosphere. Courtesy of Kevin Reardon, National Solar Observatory.

•••

5

# **Application Focus**

## The Biologist's Choice

In applications such as single molecule microscopy, super-resolution, live cell microscopy (including confocal), calcium signalling, transport/motile imaging and intracellular bioluminescence, weak, rapidly changing fluorescent signals from cells must be dynamically imaged. Andor's iXon technology offers an ideal detection solution. Ultra-sensitive detection capability in fluorescence microscopy facilitates use of lower excitation powers (thereby reducing photobleaching and phototoxicity) and lower dye concentrations.

## Single Molecule Detection

Andor's iXon EMCCD has long been the gold standard detector of the biophysics laboratory, and remains the dominant detector type, operating in a low light regime that is less suited to sCMOS cameras. The 3x accelerated 30 MHz readout speed of the iXon Ultra 888, especially combined with 'Optically Centred Crop Mode', means that dynamic single molecule processes can be better characterized. The 13  $\mu m$  pixel provides superb resolving capability at the diffraction limit.

## Spinning disk confocal

The iXon Ultra 888 is the ultimate detector to drive stunning performance from confocal spinning disk technology. Whilst affording superb confocality and low rates of phototoxicity, spinning disk experiments are inherently photon starved, by virtue of the photon rejection implicit to optical sectioning.

The superior sensitivity of the iXon Ultra detector brings these low light images to life.

#### Luminescence

The iXon Ultra can be used in either single photon counting mode (EM amplifier) or in slow scan, deep cooled CCD mode (CCD amplifier) as a highly sensitive and flexible detector in this typically long exposure, extreme low light application.


## SRRF-Stream+

Exclusive to compatible Andor cameras, SRRF-Stream leverages GPU optimization to greatly increase processing of the SRRF algorithm This makes it possible to perform super-resolution microscopy on conventional modern fluorescence microscopes in real-time!


- Real Time enhanced workflow, avoids post-processing. View in 'Live Mode'.
- ✓ Low Excitation Intensities prolonged live cell observations & accurate physiology.
- ✓ Conventional Fluorophores simple labelling, no photo-switching required.
- ✓ Live Cell Dynamics full FOV super-res images every 1-2 secs. > 10 fps using ROI.
- ✓ Cost-Effective convert conventional fluorescence microscopes to super-resolution microscopes.

Andor's new "SRRF-Stream+" provides even better performance. The original SRRF-Stream, localization was limited to 6-axes, which provided a compromise of image quality and speed for most datasets. However. for SRRF-Stream+ it has been possible to increase the axes of radiality to 24, by maximising GPU processing, enhancing image quality with minimal impact to speed. Find out more in the SRRF-Stream+ technical note.


Users of the original SRRF-Stream can upgrade to the new SRRF-Stream+ using an updater utility from their local Andor product support.

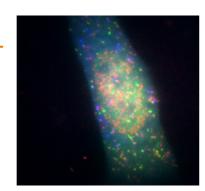

Single molecule imaging mRNA (red), during translation, and proteins, FLAG-KDM5B (green) and HA-KDM5B (blue). Courtesy of Timothy J. Stasevich, IGAF, Colorado State University.

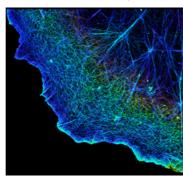

3D STORM image of actin cytoskeleton (AF647-phalloidin) in COS-7 cell. Colour codes for z-depth. Image courtesy of Xiaowei Zhuang, Harvard University.

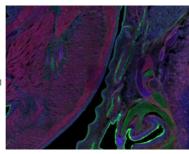

Spinning disk confocal image (right) of a mouse embryo section with F-actin (AF568-phalloidin), membrane glycoproteins (AF488-WGA), and DNA (DAPI) labelling.

## EMCCD or sCMOS?

Since the introduction of sCMOS technology by Andor, the question of the performance comparison against the more established Electron Multiplying CCD (EMCCD) has been a common one.

Being a very fast, low noise technology, sCMOS does hold some potential to offer an alternative technology across some applications and techniques, including to an extent, TIRF microscopy. Whilst the read noise of sCMOS is very low compared to CCDs, EMCCD technology holds the distinct advantage of being able to practically eliminate read noise, rendering them single photon sensitive.

After a few years of sCMOS being tested in the

staved applications for which EMCCDs were

originally purchased, such as single molecule

detection and low light spinning disk confocal

microscopy, are continuing to strongly benefit

offer a raw sensitivity that cannot be surpassed


from this ultrasensitive technology. EMCCDs

in the very low light regime

market, we are concluding that the primary light

Zyla 5.5 sCMOS 2x2 binned (13 µm)

Backilluminated EMCCD (13 µm)

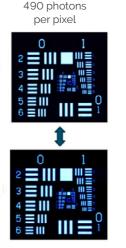

Images at a range of incident light intensity, acquired using backilluminated EMCCD iXon 888 and Zyla 5.5 sCMOS cameras (2x2 binned pixels). At low light intensities, the Signal to Noise Ratio advantage of the EMCCD is apparent (in this example the 8 photons per pixel images).



Plot of Signal to Noise Ratio versus Incident Photon Intensity, comparing a back-illuminated EMCCD iXon 888 (13  $\mu$ m pixel size) to 2x2 binned Zyla sCMOS cameras (13  $\mu$ m pixel size after binning)

6 7

## **Technical Specifications**

System Specifications •2

	Ultra	888		Ultra 897				
Sensor QE options	#BV: Back Illuminated, standard AR coated UVB: Back Illuminated, standard AR with additional lumogen coating EXF: Back illuminated, dual AR coated with fringe suppression #BB: Back-illuminated, blue optimized AR coated							
Fringe Suppression		Availab	ole on EXF sensor optio	n				
Active pixels	1024)	< 1024		512 x 512				
Pixel size	13 x 1	13 μm		16 x 16 μm				
Image area	13.3 x 13.3 mm w	ith 100% fill factor	8.2 x 8.2	mm with 100% fill factor				
Pixel Readout Rate Minimum temperature, air cooled, ambient 20°C Chiller liquid cooling, coolant @ 10°C, >0.75l/min	10 MHz -80°C -95°C	30 MHz <sup>•3</sup> -60°C -75°C	10 MHz -80°C -100°C	17 MHz -80°C -100°C				
Thermostatic Precision			± 0.01°C					
Triggering	Inte	ernal, External, Externa	l Start, External Exposu	ıre, Software Trigger				
System window type	UVB, EXF: UV-grad	de fused silica, Broadb	and Vacuum Ultraviole	nfrared, 0.5 degree wedge et-Near Infrared, 0.5 degree wedge Near Infrared, 0.5 degree wedge				
Blemish specification	Grade 1	sensor from supplier.	Camera blemishes as c	lefined by Andor Grade A				
Digitization		16-bit (at all speeds)						
PC Interface	USB 3.0 <sup>•12</sup> USB 2.0							
Lens Mount		C-mount						
Direct Data Access		Cam	era Link 3-tap output					

## Advanced Performance Specifications •2

			Uli	ra 88	8				Uli	tra 897			
Dark current and background events <sup>e4,5</sup>													
Dark current (e-/pixel/sec) @ -80°C Dark current (e-/pixel/sec) @ max cooling	0.00025 0.00011				0.00030 0.00015								
Spurious background (events/pix) @ 1000x gain / -85°C	0.005			0.0018									
Active area pixel well depth			80	,000 €	<b>&gt;</b> -				180	0,000 e			
Gain register pixel well depth <sup>6,7</sup>	730,000 e <sup>-</sup>				800,000 e <sup>-</sup>								
Pixel readout rates	EM Amplifier: 30, 20, 10 & 1 MHz Conventional Amplifier: 1 & 0.1 MHz			EM Amplifier: 17, 10, 5 & 1 MHz Conventional Amplifier: 3, 1 & 0.08 MHz									
Read noise (e-) <sup>e7</sup>			CCD olifier			entional aplifier					onventional Amplifier		
MHz	30	20	10	1	1	0.1	17	10	5	1	3	1	0.08
Without Electron Multiplication With Electron Multiplication	130	80	40	12	6	3.5	89	65	37	15	9.6	5.3	2.7
with Electron Multiplication	< 1	< 1	< 1	< 1	-	-	< 1	< 1	< 1	< 1	-	-	-
Linear absolute Electron Multiplier gain		1 - 10	000 tin	nes vid	a Real(	Gain™ (cali	bration	stable	at all co	ooling te	mperc	itures)	)
Linearity*8						Bette	er than 9	99.9%					
Vertical clock speed	0.6 to 4.33 µs (user selectable)					0.3	3 to 3.33	3 µs (use	r sele	ctable	)		
Timestamp accuracy	10 ns												
NEW SRRF-Stream* mode						(	Optiono	al					

## iXon Ultra 888 Frame Rates

## Standard Mode \*3,9

Binning	1024 x 1024	512 x 512	256 x 256	128 x 128	1024 x 100	1024 x 32	1024 x 1
1 x 1	26	50	95	171	220	498	1163
2 x 2	50	94	170	285	368	699	-
4 x 4	92	167	281	426	552	870	-

## Crop Mode - Optically Centred frame rates in brackets • 3.9

Binning	512 x 512	256 x 256	128 x 128	64 x 64	1024 x 100	1024 x 32	1024 x 1
1 x 1	93 (78)	190 (251)	670 (697)	2053 (1319)	259	778	9690
2 x 2	170 (143)	350 (426)	1150 (1019)	3123 (1646)	492	1416	-
4 x 4	291 (245)	601 (653)	1772 (1504)	4109 (1857)	887	2370	-

## iXon Ultra 897 Frame Rates

## Standard Mode •10


Binning	512 x 512	256 x 256	128 x 128	64 x 64	512 x 100	512 x 32	512 x 1
1 x 1	56	110	212	398	267	708	2,881
2 x 2	109	210	394	699	486	1,141	-
4 x 4	206	385	682	1,109	820	1,615	-


## Crop Mode - Optically Centred frame rates in brackets<sup>10</sup>

Binning	256 x 256	128 x 128	64 x 64	32 x 32	512 x 100	512 x 32	512 x 1
1 x 1	111 (174)	595 (569)	1,433 (1,490)	3,533 (3,021)	282	857	11,074
2 x 2	215 (329)	1,094 (1,013)	2,481 (2,325)	5,555 (4,048)	541	1,607	-
4 x 4	405 (593)	1,883 (1,661)	3,906 (3,236)	7,751 (4,878)	1,005	2,865	-

## Quantum Efficiency (QE) Curves \*11

# Blue-optimized sensor option for quantum and BEC studies

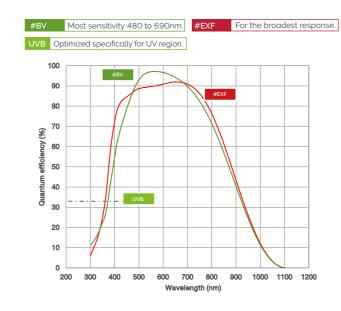

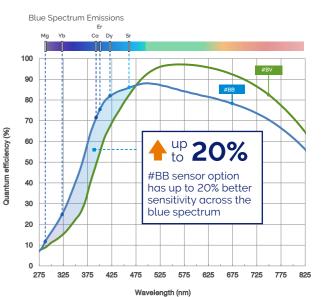

## Creating The Optimum Product for You

## Step 1.

#### Choose the camera type



Description	Code
iXon Ultra 888: 1024 x 1024 EMCCD, max. 30 MHz, with USB 3.0	DU-888U3-CSO-
iXon Ultra 897: 512 x 512 EMCCD, max. 17 MHz, with USB 2.0	DU-897U-CSO-
Add sensor type code to camera code (see step 2)	

Type

NOTE: If SRRF-Stream\* real time super-resolution functionality is required with your iXon Ultra, please order as an 'accessory' in step 4 below.

### Step 2.

#### Choose the sensor type option



Type

Description	Code
Back-illuminated, standard AR coated	#BV
Back Illuminated, standard AR with additional lumogen coating	UVB

Description						
Back-illuminated, EX2 dual AR coated with fringe suppression	EXF					
NEW Back-illuminated, blue optimized AR coated	#BB					

Add sensor type code to camera code (step 1)

#### Step 3.

### Select an alternative camera window (optional)



The standard window has been selected to satisfy most applications. However, other options are available. The alternative camera window code must be specified at time of ordering.

To view and select other window options please refer to the 'Camera Windows Supplementary Specification Sheet' which gives the transmission characteristics, product codes and procedure for entering the order. Further detailed information on the windows is in the Technical note - 'Camera Windows: Optimizing for Different Spectral Regions'

## Step 4.

#### Select the required accessories

	Description	Order Code	Description	Order Code
Accessories	SRRF-Stream+ real time super-resolution functionality, compatible with iXon Ultra and iXon Life EMCCD platforms. Camera must be connected to acquisition PC workstation containing an Nvidia GPU card (compute capability v3.0, or above, and 4GB or areater on-board GPU RAM).	SRRF- STREAM- IXON	Re-circulator for enhanced cooling performance	XW-RECR
			Oasis 160 Ultra compact chiller unit (tubing to be ordered separately)	ACC-XW- CHIL-160
	SRRF-Stream Dell Workstation (English), pre-installed with a recommended and tested GPU card, alongside SRRF-Stream enabled MicroManager and Andor SDK2 with SRRF-Stream.	WKST-SRRF- 9ZY	6 mm tubing options for ACC-XW-CHIL-160 (2x2.5 m or 2x5m lengths)	ACC-6MM- TUBING-2X2.5/ ACC-6MM- TUBING-2X5M
	Monitor (optional) - Dell UltraSharp U3417W - 34.14" Curved LED	FUS-MNTR- 34W	C-mount to Nikon F-mount adapter	OA-CNAF
			C-mount to Olympus adapter	OA-COFM
	Dell UltraSharp UP3017 - 30" with PremierColor	FUS- MNTR-30	C-mount to T-mount adapter	OA-CTOT
	OptoMask accessory, used to mask unwanted sensor area during Crop Mode acquisition (refer to OptoMask Specification Sheet for further information).	OPTMSK-L/ OPTMSK- OC-L/ OPTMSK- OC-S	15 m Active USB 3.0 connector cable (power supply not required) Icron for Ultra 888	ACC-ASE-06887
			50 m Fibre Optic USB 3.0 extender solution inc. power supply (Adnaco) for Ultra 888	ACC-ASE-08762
			100 m Fibre Optic USB 3.0 extender solution inc. power supply (Adnaco) for Ultra 888	ACC-ASE-07860

### Step 5.

#### Select the required software

The iXon Ultra series requires one of the following software options:

Solis Imaging: A 32-bit and fully 64-bit enabled application for Windows (8, 8.1 and 10) offering rich functionality for data acquisition and processing. AndorBasic provides macro language control of data acquisition, processing, display



Andor SDK: A software development kit that allows you to control the Andor range of cameras from your own application. Available as 32 and 64-bit libraries for Windows (8, 8.1 and 10), compatible with C/C++, C#, Delphi, VB.NET, LabVIEW and Matlab. Linux SDK compatible with C/C++.

Andor iQ A comprehensive multi-dimensional imaging software package. Offers tight synchronization of EMCCD with a comprehensive range of microscopy hardware, along with comprehensive rendering and analysis functionality. Modular architecture for best price/performance package on the market.

Third party software compatibility, drivers are available for a variety of third party imaging packages

For SRRF-Stream the iXon must be operated either through MicroManager (Open Imaging) open source microscopy software platform, or through the Andor SDK, if SRRF-Stream functionality is to be accessed.

# **Product Drawings**


Dimensions in mm [inches]


#### iXon Ultra 888


# 139.6 [5.496] MOUNTING POINTS 1/4 - 20 UNC X 12.7 [0.500] DEEP 6 PLCS 46.5[1.831] 46.5[1.831]


## 139.6[5.496] 120.0[4.724] 46.5 [1.831] 46.5 [1.831] 38.1[1.500] 38.1[1.500]


iXon Ultra 897

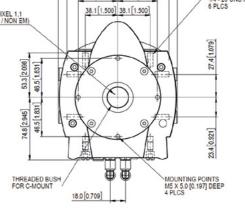

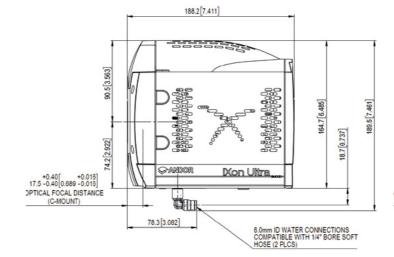

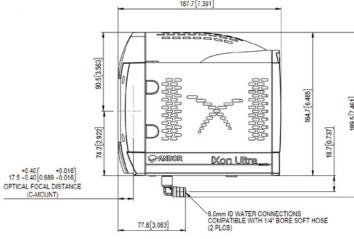

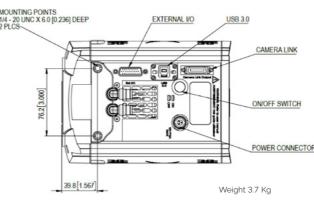

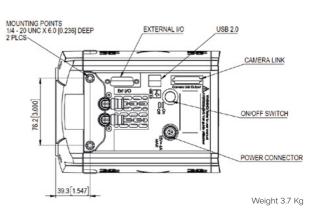

## **Ultra 888 Power Requirements**

- Power Input: +12 VDC ± 5% @ 8 A
- Power Consumption: 96 W max
- Ripple and noise: 120 mV max (peak-peak 0 20 MHz)
- External Power Supply: 100 240 VAC 50/60 Hz

#### **Ultra 897 Power Requirements**

- Power Input: +12 VDC ± 5% @ 6 A
- Power Consumption: 72 W max
- Ripple and noise: 120 mV max (peak-peak 0 20 MHz)
- External Power Supply: 100 240 VAC 50/60 Hz

Logic: Connector type: 26 way D Type with 8 programmable digital inputs or outputs for control and sensing of up to 8 external devices, Minimum cable clearance required: 90 mm, Weight: 3.7 kg [8 lb 3 oz] approx.

10 11



# Order Today

Need more information? At Andor we are committed to finding the correct solution for you. With a dedicated team of technical advisors, we are able to offer you one-to-one guidance and technical support on all Andor products. For a full listing of our local sales offices, please see: andor.oxinst.com/contact

### Our regional headquarters are:

## **Europe**

Belfast, Northern Ireland Phone +44 (28) 9023 7126 Fax +44 (28) 9031 0792

#### **North America**

Concord, MA, USA Phone +1 (860) 290 9211 Fax +1 (860) 290 9566

#### Japan

Tokyo Phone +81 (3) 6732 8968 Fax +81 (3) 6732 8939

#### China

Beijing Phone +86 (10) 5884 7900 Fax +86 (10) 5884 7901

## **Items shipped with your iXon Ultra 888:** 1x Andor ACZ-03463: 2m Multi I/O timing

cable, offering Fire, External Trigger, Shutter and Arm

1x 3m USB 3.0 cable Type A to Type B PCIe USB 3.0 Card Adapter (2-Port)

1x Power supply unit with mains cable

1x Quick Start guide

1x Electronic copy of user manuals 1x SRRF-Stream Quick Start guide (if applicable)

1x Individual system performance booklet

Items shipped with your iXon Ultra 897: 1x Andor ACZ-03463: 2 m Multi I/O timing cable, offering Fire, External Trigger, Shutter

and Arm 1x 3m USB 2.0 cable Type A to Type B

1x Power supply unit with mains cable 1x Quick Start guide

1x Electronic copy of user manuals
1x SRRF-Stream Quick Start guide (if applicable)

1x Individual system performance booklet

#### **Recommended Computer Requirements:**

- 3.0 GHz single core or 2.6 GHz multi core processor
- 100 MB free disc space to install software (at least 1 GB recommended for data spooling)
- USB 3.0 Super Speed Host Controller capable of a sustained rate of 60MB/s for iXon Ultra 888
- USB 2.0 High Speed Host Controller capable of sustained rate of 40MB/s for iXon Ultra 897
- Solid-state drive (SSD) capable of a minimum sustained write speed of 100MB/S for spooling data
- Windows (8.1 and 10) or Linux
- SRRF-Stream+ If selected, the PC requires a Nvidia GPU card. See page 10 for further

## Footnotes: Specifications are subject to change without notice

- Assembled in a state-of-the-art cleanroom facility, Andor's
- vacuum seal (no o-rings), with a stringent protocol to minimize outgassing, including use of proprietary materials Figures are typical unless otherwise stated.
- At 30 MHz overclocked pixel readout rate, thermal dissipation from the sensor is higher since a greater proportion of time is spent vertical shifting, and it is necessary to set a higher sensor cooling temperature at this rate. Furthermore, stable cooling performance will depend on other variables such as vertical clock speed. Region of Interest size (Standard or Crop Mode) and ambient temp. As such, user testing is advised to determine the stable sensor cooling temperature for
- selected conditions. Status of temperature stability is apparent through the acquisition software. The dark current measurement is averaged over the sensor area excluding any regions of blemishes
- The dark current measurement is averaged over the sensor area excluding any regions of blemishes. Using Electron Multiplication the iXon is capable of detecting single photons, therefore the true camera detection limit is set by the number of dark' background events. These events consist of both residual thermally generated electrons and Clock Induced Charge (CiC) electrons (also referred to as Spurious Noise), each appearing as random single spikes above the read noise floor. A thresholding scheme is employed to count these single electron events and is quoted as a probability of an event per pixel. Acquisition conditions are full resolution and max frame rate (30 MHz readout; frametransfer mode: 1.1 µs vertical clock speed; x 1000 EM gain; 10 ms exposure; -95°C).

  The EM register on CCD201 sensors has a linear response up to -400,000 electrons and a full well depth of -730,000 electrons.
- Readout noise is for the entire system. It is a combination of sensor readout noise and A/D noise. Measurement is for Single Pixel readout with the sensor at a temperature of -75°C and minimum exposure time under dark conditions. Under Electron Multiplying conditions, the effective system readout noise is reduced to sub 1 e<sup>-</sup> levels. Linearity is measured from a plot of counts vs. exposure time under constant photon flux up to the saturation point of the system, at 10 MHz readout speed.

  All measurements are made at 30 MHz pixel readout speed with 0.6 µs vertical clock speed. It also assumes internal trigger made of postpration. Strangard and Crop Mede frame rates shown are for "Corner Tethorical" Pole with "Obtically trigger made of postpration."
- trigger mode of operation. Standard and Crop Mode frame rates shown are for 'Corner Tethered' ROIs, with 'Optically Centred' ROI frame rates shown within brackets.
- 10.All measurements are made at 17 MHz pixel readout speed with 0.3 µs vertical clock speed. It also assumes internal trigger mode of operation. Standard and Crop Mode frame rates shown are for 'Corner Tethered' ROIs, with 'Optically
- Centred ROI frame rates shown within brackets.

  11. Quantum efficiency of the sensor at 25°C, as supplied by the sensor manufacturer.

  12. IXon Ultra 888 should work with any modern USB 3.0 enabled PC/laptop, as every USB 3.0 port should have its own host controller. IXon Ultra 888 also ships with a USB 3.0 PCI card as a means to add a USB 3.0 port to an older PC, or as a diagnostic aid to interoperability iss

#### **Operating & Storage Conditions**

- Operating Temperature: -20°C to 30°C ambient
- Relative Humidity: < 70% (noncondensing)
- Storage Temperature: -25°C to 50°C

#### **Power Requirements**

• Please refer to page 11

















