

Configure Sona or Marana with µManager

ANDOR

Covers Sona and Marana Models 4.2B-11, 4.2B-6, 2.0B-11, USB 3.0 and CoaXPress variants

andor.oxinst.com

Sona or Marana with Micro-Manager

Part 1

Install the latest Micro-Manager build



Install Micro-manager 2.0 beta nightly from here: https://micro-manager.org/wiki/ Version_2.0#Downloads

Version 2.0

Micro-Manager 2.0 is a major step forwards compared to the 1.4 series that preceded it. The user interface (UI) has been overhauled, especially with respect to the image display window. The API has also been completely reworked, allowing for more powerful scripts and plugins. This was all enabled by a new internal data model for working with images and image metadata. This page serves as a "hub" for various other pages related to Version 2.0. Contents 1 Downloads
2 Contributing to MM2.0
3 API
4 API Events
5 Other Documentation

Downloads


Beta nightly builds: Windows | Mac OS X

2.0.0-beta2 (obsolete) Windows 32-bit | Windows 64-bit | Mac OS X (Note: there is no 2.0.0-beta1.) Alpha nightly builds (obsolete): Windows | Mac OS X) Current source code on GitHub: mm2 branch

Contributing to MM2.0

How to debug and develop MM2.0

Follow the setup wizard steps:

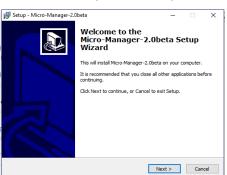

Index of /~MM/nightlyBuilds/2.0.0-beta/Windows

	Name	Last modified	Size	Description
	Parent Directory			
	Parent Directory		-	
	MMSetup 64bit 2.0.0-beta3 20190228.exe	28-Feb-2019 23:01	70M	
101 10	MMSetup 32bit 2.0.0-beta3 20190228.exe	28-Feb-2019 23:01	68M	
10 10 10	MMSetup 64bit 2.0.0-beta3 20190227.exe	27-Feb-2019 23:02	701	
	MMSetup 32bit 2.0.0-beta3 20190227.exe	27-Feb-2019 23:02	68M	
191 010	MMSetup 64bit 2.0.0-beta3 20190226.exe	26-Feb-2019 23:00	70M	
181 181	MMSetup 32bit 2.0.0-beta3 20190226.exe	26-Feb-2019 23:00	68M	
10	MMSetup 64bit 2.0.0-beta3 20190225.exe		70M	
	rmsetup 04010 2.0.0-Deta3 20190225.exe	25-rep-2019 23:00	761	
10- 10- 10-	MMSetup 32bit 2.0.0-beta3 20190225.exe	25-Feb-2019 23:00	68M	

- 1. The latest build will be at the top of the list. Example here shows build for 28-Feb-2019
- 2. Admin rights are required to run the executable file
- 3. The Micro-Manager install file directory path is needed for Part 2

Sona or Marana with Micro-Manager

Install the Andor Driver Pack 3 Download the following exe file Andor Driver Pack - min version 3.15 (sCMOS) at andor.oxinst. com/downloads Install the Andor Driver Pack 3 into to the Micro-Manager install directory (from Part 1) Extract Compressed (Zipped) Folders Select a Destination and Extract Files Files will be extracted to this folder C:\Program Files\Micro-Manager-2.0beta Browse... Show extracted files when complete Extract Cancel Additional files will now be present in the Micro-Manager install directory, an example of some of these additional files is shown below: 🗟 atcore.dll 05/03/2019 13:50 atdevchamcam.dll 05/03/2019 13:50

Notes

1. The install directory should be set during the micro-manager install, in this example default was: C:/Program Files/Micro-Manager-2.0beta

2. The Andor Driver Pack will be updated regularly and the version number of the file may differ from that given above. The most recent release can be found at <u>andor.oxinst.com/downloads/</u>



Help and Support

Please see our range of technical documentation available at: <u>andor.oxinst.com/downloads</u> please create an account to request access to software files.

If you require further help please contact our customer support team at andor.oxinst.com/support

Further help on third party software compatibility can be found in our third party software compability matrix.

.